

# MilliKey™ 24mm Button Caps

MilliKey MH models can optionally use different interchangeable solid color and clear-top button caps, which can be purchased in packs including 8 to 12 button caps and a removal pin.



## Color Cap Pack (24mm)

CP24-CMIX-12

Includes 12 solid color caps (2 red, 2 yellow, 2 green, 2 blue, 2 black, and 2 pink) and one cap removal pin.



## Clear Cap Pack (24mm)

CP24-CLKW-8

Includes 8 clear caps, four with white and four with black inner tops, and one cap removal pin.




For further information or to place an order, please visit our website or email [sales@labhackers.com](mailto:sales@labhackers.com).




**LabHackers Inc.**

[www.labhackers.com](http://www.labhackers.com)


# Changing MilliKey™ Button Caps

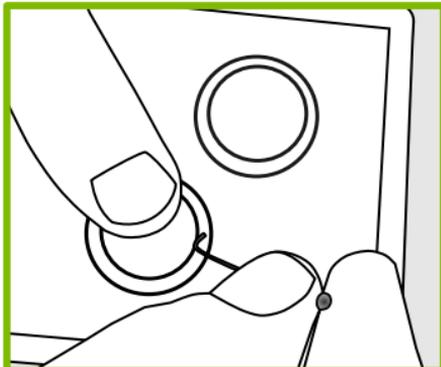


**1** Insert removal pin between button cap and outer button ring.

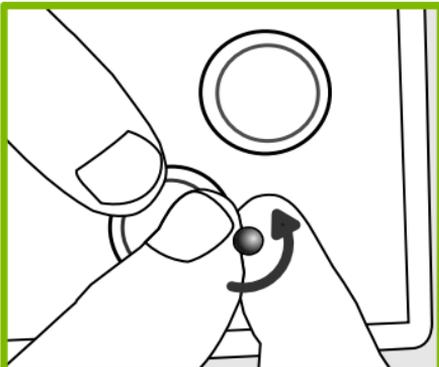

**2** Push pin to button bottom and rotate ~ 45° towards button cap.

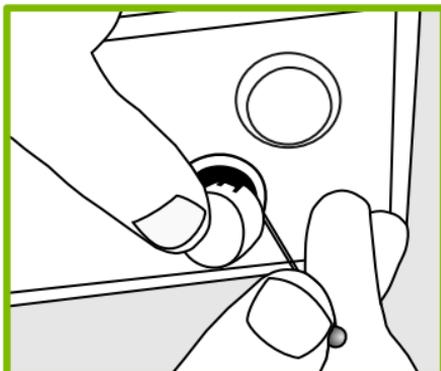

**3** Place finger over cap, pull up on pin to lift cap out of the box.

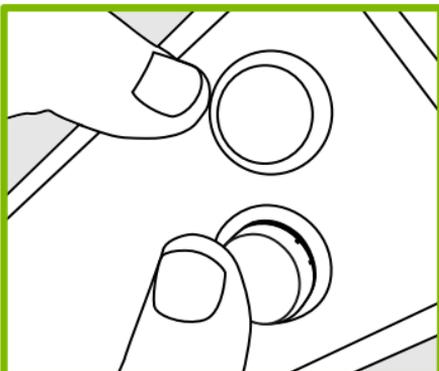

**4** Insert new cap into button hole and rotate until cap is in place.

Please email [support@labhackers.com](mailto:support@labhackers.com) with any questions.



**LabHackers Inc.**

[www.labhackers.com](http://www.labhackers.com)