

Quickstart Guide

MilliKey™ MH-5 Response Box



Default MilliKey MH-5 Button to Keyboard Key Mappings

Thank you for choosing MilliKey.

The MilliKey MH-5 is a 1000 Hz USB Response Box that generates standard keyboard events when buttons are pressed.

Initial Set Up

■ Keyboard

The MilliKey is automatically detected as a 1000 Hz USB HID keyboard when connected to a Windows, macOS, or Linux computer. Start collecting standard keyboard data, it is that simple.

■ Serial

MilliKey response boxes also optionally utilize a 1000 Hz USB Serial interface, which is plug-and-play on Windows 10, macOS, and Linux. If using Windows 7 please install the required USB Serial driver from:

www.labhackers.com/downloads.html



LabHackers Inc.

www.labhackers.com

Quickstart Guide cont'd


MilliKey™ MH-5 Response Box

Modifying Settings

The LabHackers Device Manager application makes it easy to view and modify MilliKey settings such as button to keyboard key mappings.

Download Device Manager for Windows from:

www.labhackers.com/downloads.html

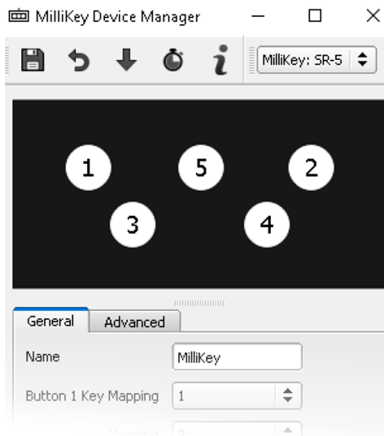

Further Information

For any questions about the MilliKey MH-5, please contact:

Sales: sales@labhackers.com

Support: support@labhackers.com

For more information about other MilliKey models and options, please visit www.labhackers.com.



LabHackers Inc.

www.labhackers.com